



Product Usage

Product: **Chicken Embryo Extract**
Powder

Catalog #: 100-163P

Product Description

Gemini's chicken embryo extract is prepared from 9-12 day old eggs. It is processed as a 50% extract in phosphate-buffered saline that is supplemented with antibiotic:antimycotic. The liquid extract is sterile filtered to 0.2 μm and lyophilized in 10mL aliquots. The sterility of this lyophilized powder cannot be guaranteed and should be sterile filtered once more after reconstitution.

Storage

Recommended storage is -20°C or below. Lyophilized product may be stored at $2-8^{\circ}\text{C}$.

Directions for Use

Reconstitute lyophilized product to 10ml using sterile water. Agitate to mix or use an orbital shaker set to 150 rpm and mix for 30 minutes.

It is advised that lyophilized product is filtered to 0.2 μm after reconstitution to maintain sterility. For best filtering results, use a pre-filter then 0.45 μm and finish with 0.2 μm filter.

For *in vitro* diagnostic laboratory use or further manufacturing only. Not for human therapeutic use.